SNIFFY, O RATO VIRTUAL

CONDICIONAMENTO OPERANTE

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Sniffy

the Virtual Rat, "Pro Version"

Welcome to our virtual laboratory for studying the learning and behavior of Sniffy, the Virtual Rat,

PRO VERSION

The **Overview** illustrates the main features.

The **Projects** illustrate how one might use Sniffy Pro to teach Operant and Classical conditioning and Discrimination and Generalization.

Please choose one of the options below.

Overview

Operant Conditioning

Project 2

Project 3

Exit

Condicionamento Operante

About


Sniffy

the Virtual Rat, "Pro Version"

Project 1:

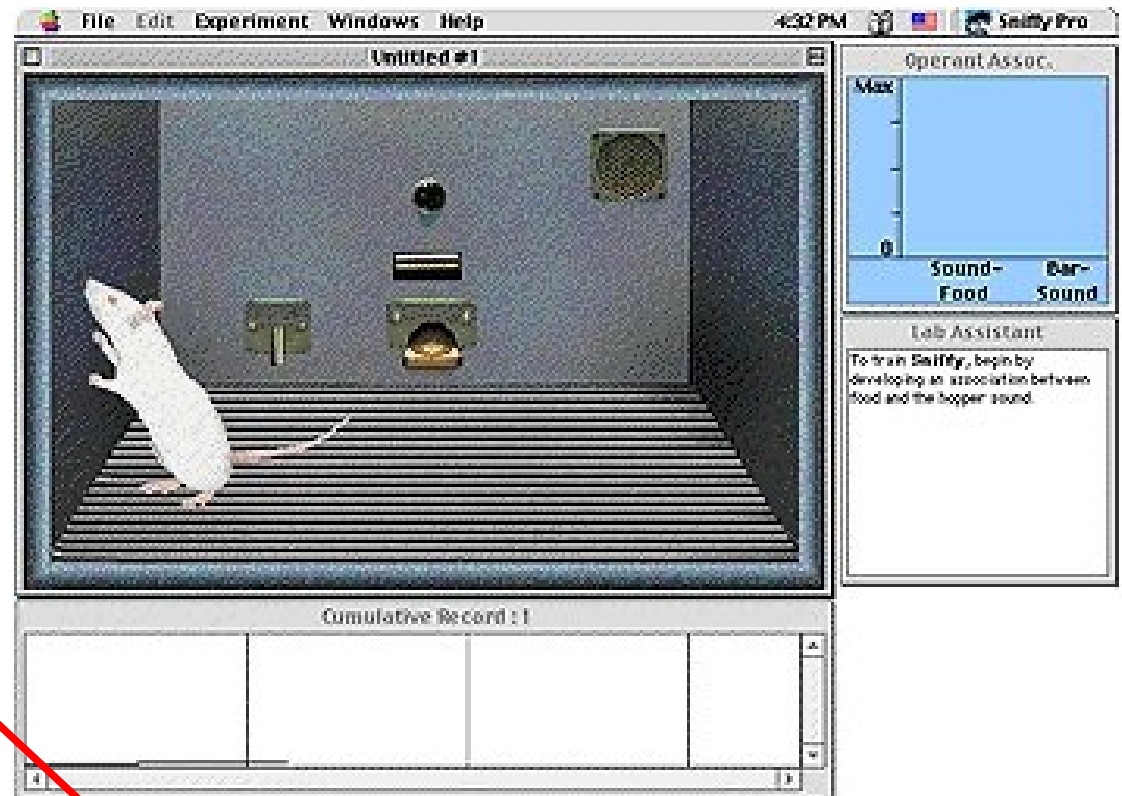Basic Operant Phenomena:
Magazine Training, Shaping,
Extinction and Spontaneous
Recovery.

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Stage 1: Magazine Training. When you start the program you will see Sniffy exploring the chamber. To train Sniffy to press the bar, you must first present food pellets so that Sniffy learns to associate the sound of the magazine with the availability of food.

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Wait until Sniffy closely approaches the food hopper. Then deliver a food pellet either by pressing the space bar on your computer keyboard or by clicking your mouse button while pointing at the bar.

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
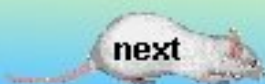
You may want to give Sniffy several pellets in rapid succession before he wanders away from the hopper. After Sniffy has received several pellets, you can let him wander away a short distance before giving him the next pellet.

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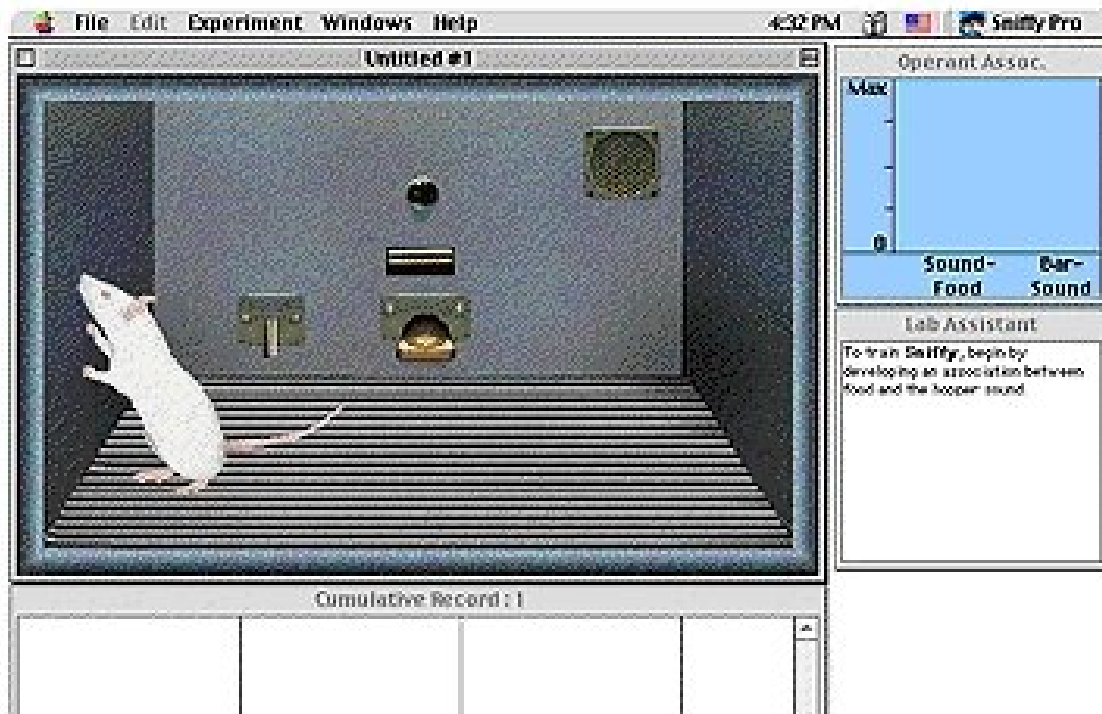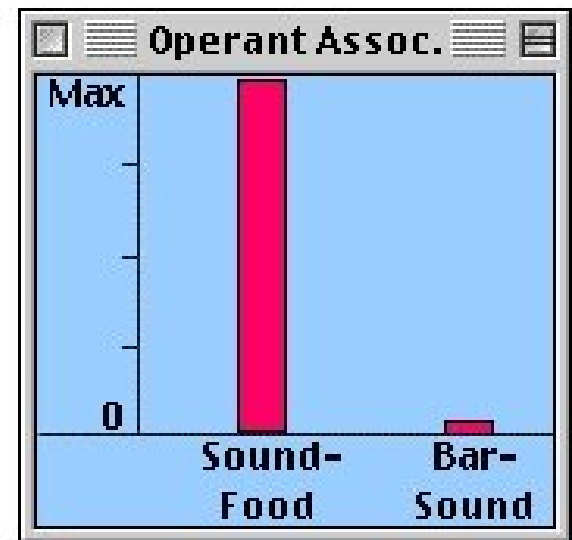
Lab Assistant

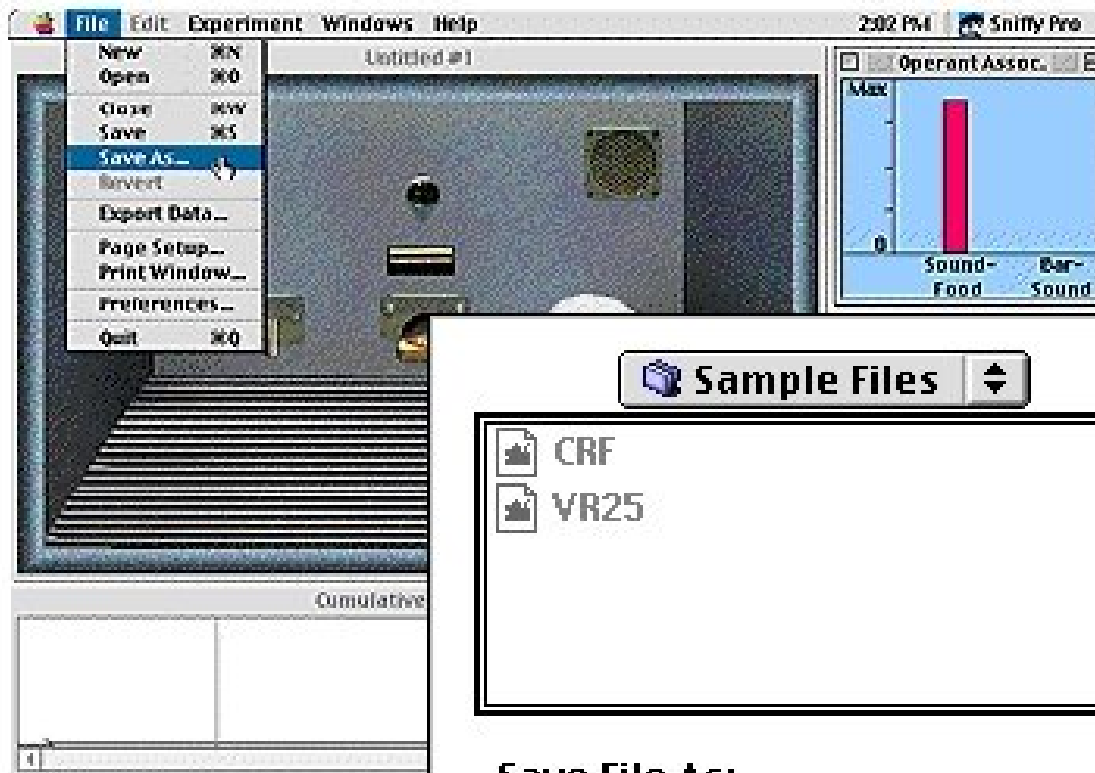
Sniffy appears to have developed an association between the sound of the hopper and the food. You can now use the sound as a reinforcer to increase **Sniffy's** rearing behaviours. Once you have trained **Sniffy** to rear often, begin to refine the behaviour by only reinforcing rearing near the hopper. **Sniffy** should learn to bar press.

Sniffy is magazine trained when the vertical bar for Sound-Food in the Operant Associations mind window reaches the maximum. At that point, no matter where Sniffy is, he will go get the food pellet whenever he hears the sound of the magazine (the pellet dispenser).

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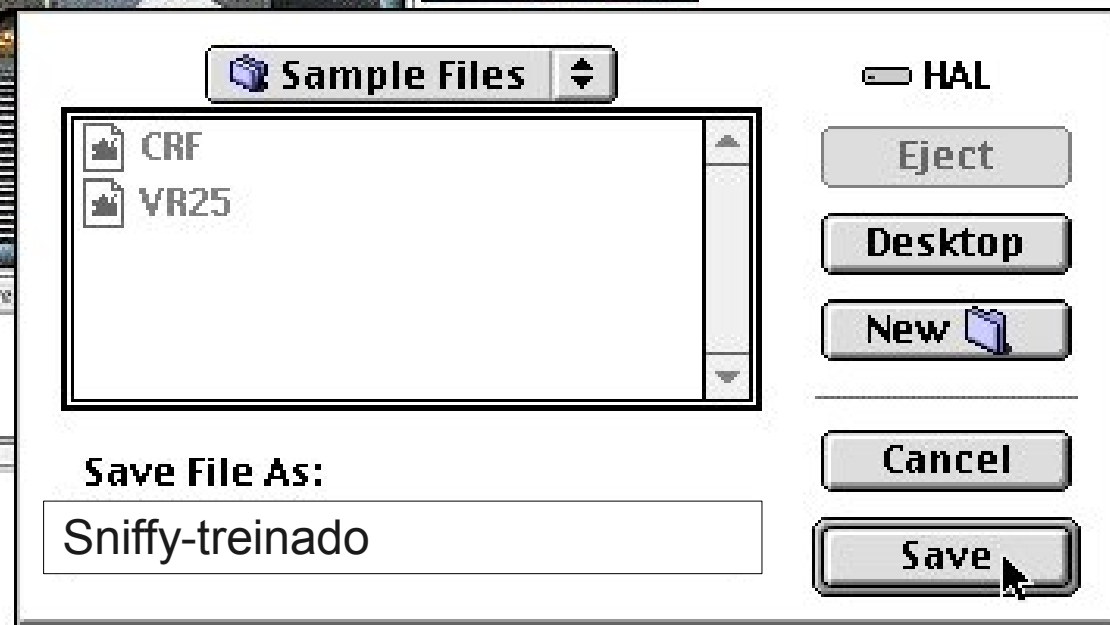
You should save a file after each stage of training. Select the Save As command from the File menu, give the file a name (e.g. MagTrain), and save it in the folder where you want it (e.g. Sniffy Pro folder).

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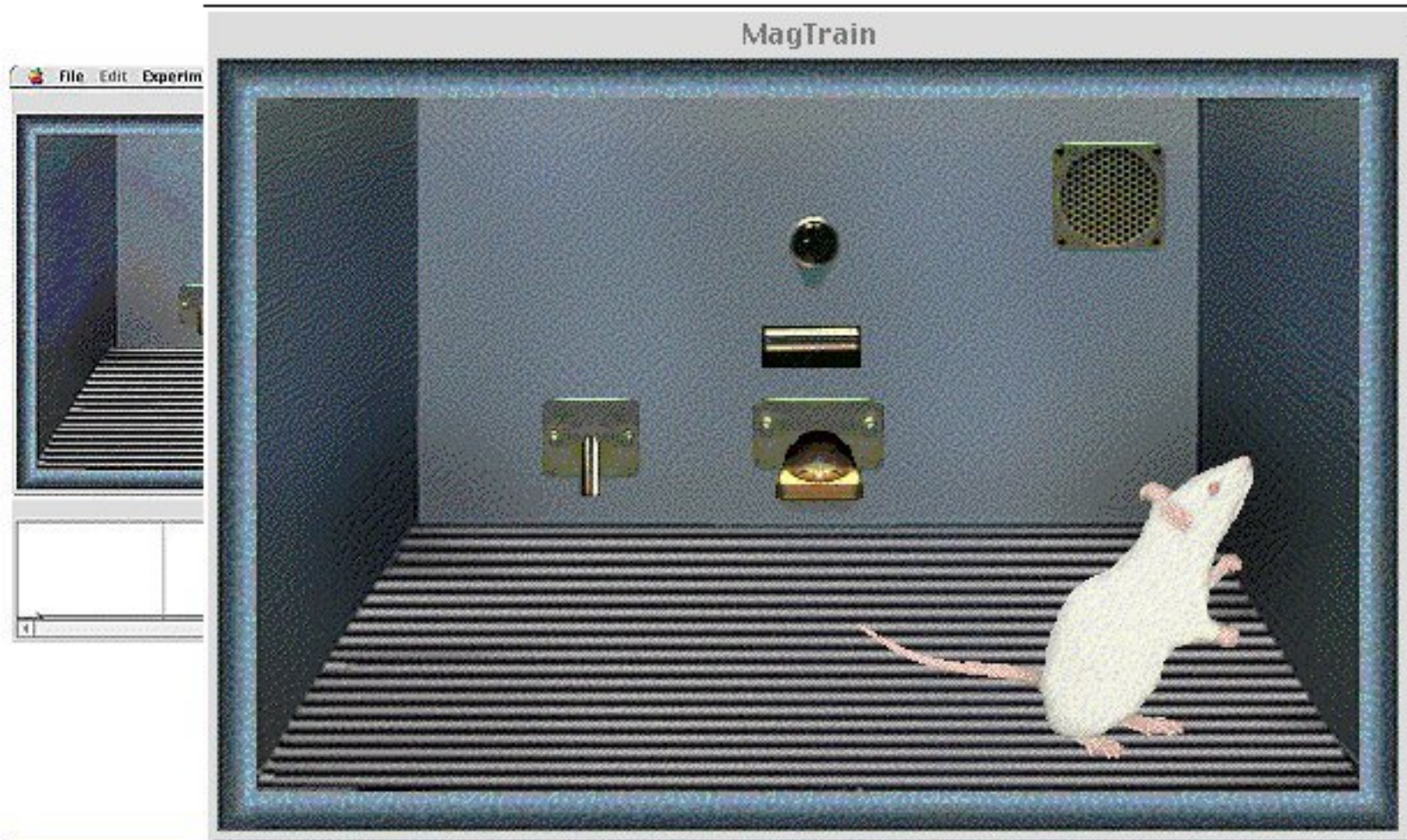
Shaping Sniffy to press the bar. To begin shaping, reinforce Sniffy when he rears up anywhere in the operant chamber. Timing is critical. Click on the bar or press the spacebar to present food **while** Sniffy is rearing up.

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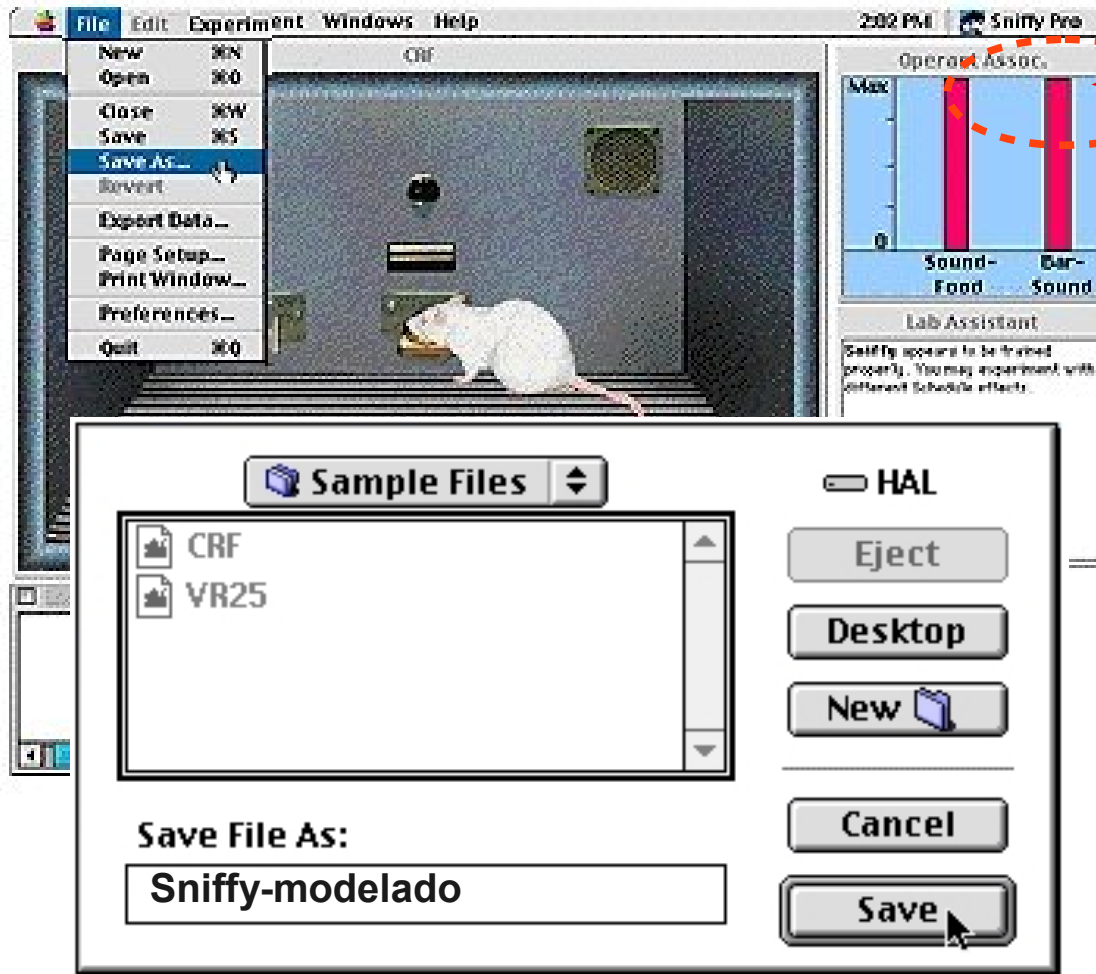
Once rearing has become more common; reinforce Sniffy only when he rears up against the back wall where the bar is located. Gradually require Sniffy to rear up closer and closer to the bar.

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When Sniffy is fully trained, select the Save As... command from the File menu and name your file to preserve your trained Sniffy for future use.

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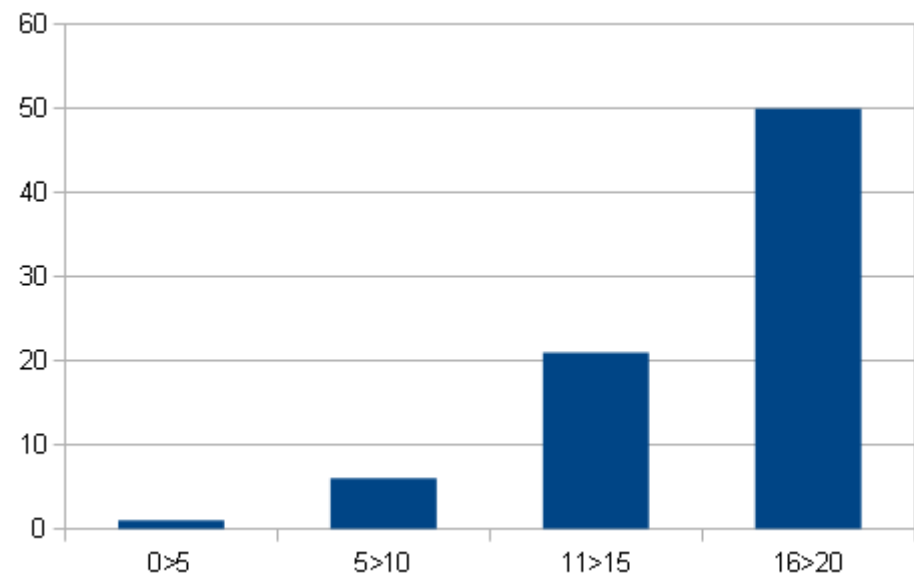
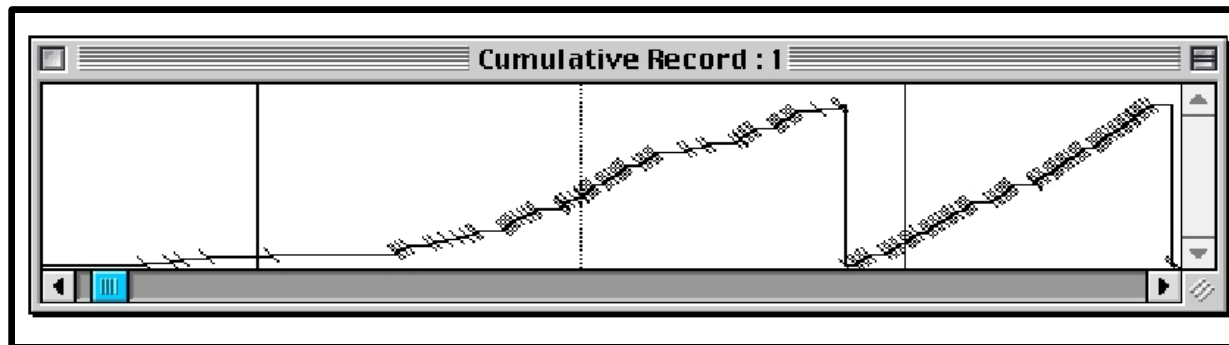
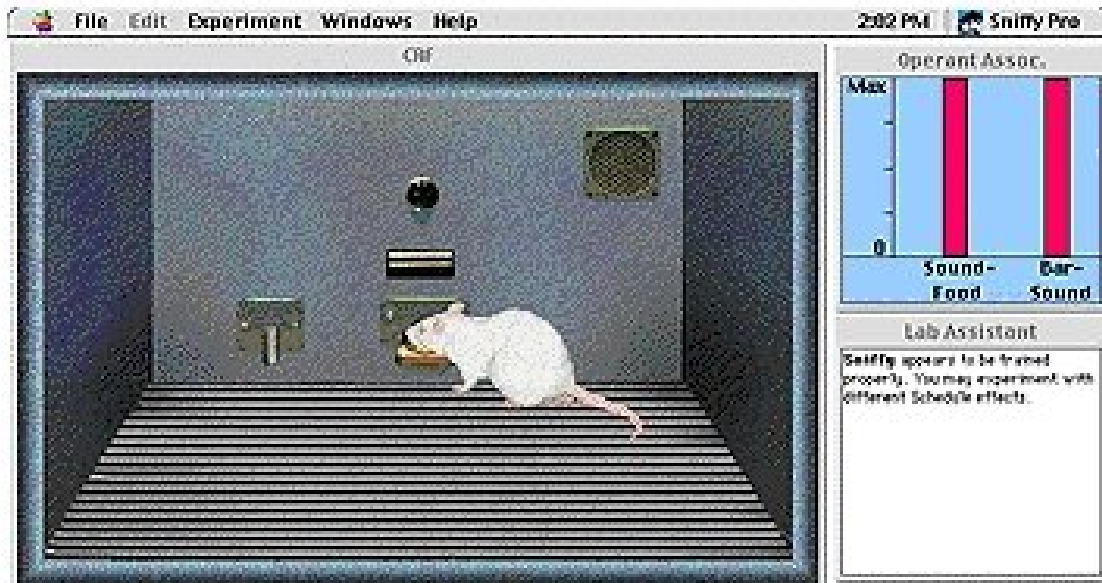
→ Menu → Windows → Response Measures → Cumulative Record

The Cumulative Record is normally shown under the Operant Chamber window. If it has been closed or it is not visible, you can make it visible by selecting it from the Windows menu.

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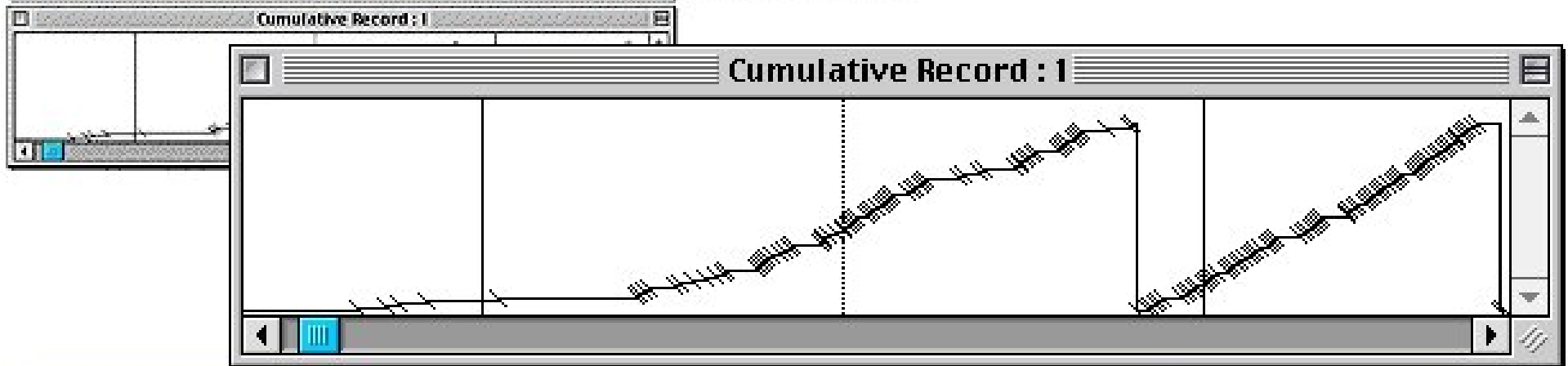
Changes in the bar-pressing rate are measured in a graph called the **Cumulative Record**. The horizontal dimension of the graph represents time, and the vertical dimension represents bar presses.

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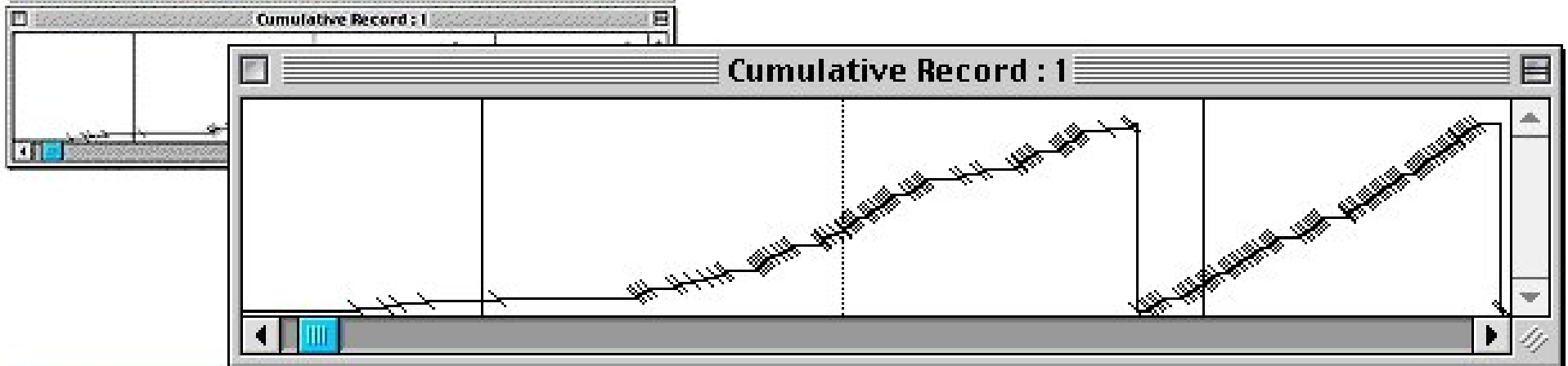
As time elapses a line is drawn horizontally across the page. Each time Sniffy presses the bar the line moves up a notch and continues along horizontally. Reinforced responses are marked by short, oblique lines drawn through the record.

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
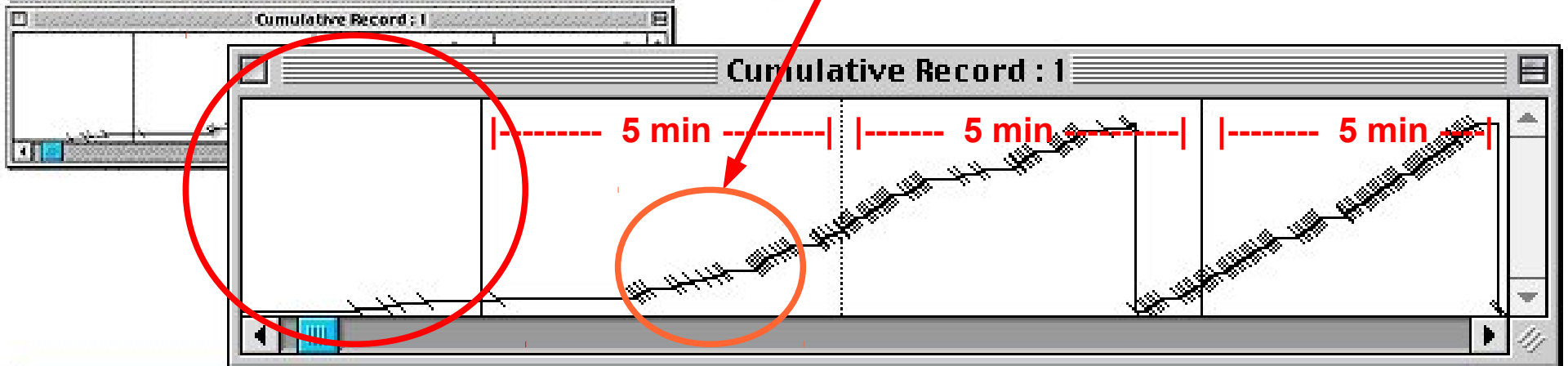
On the left, the horizontal line shows that Sniffy has not yet learned to press the bar. Sniffy begins to press the bar slowly at first and then more and more frequently. The dark vertical lines are pen resets from the top to the bottom. The thin, solid and dotted vertical lines are spaced at 5-minute intervals.

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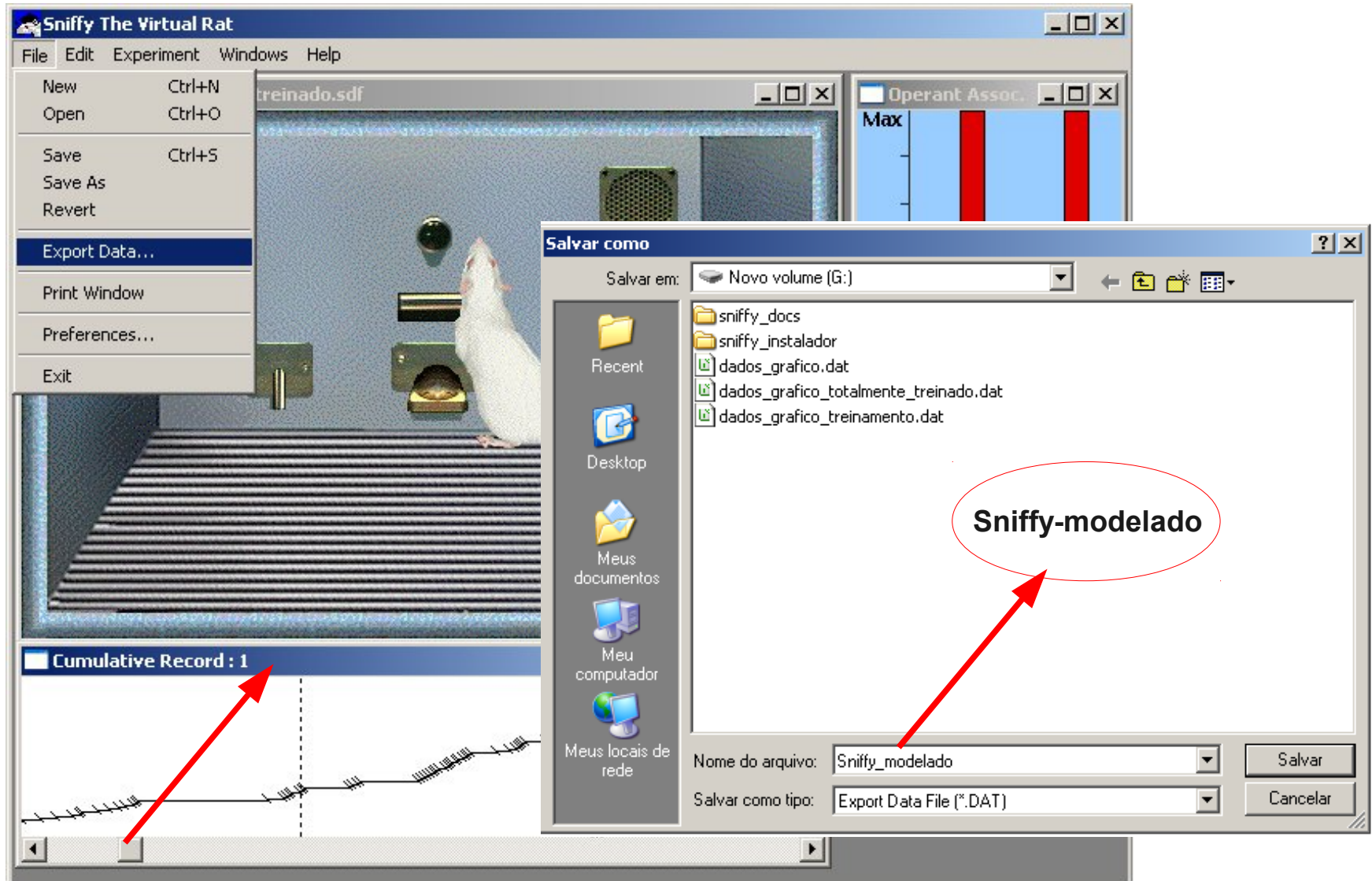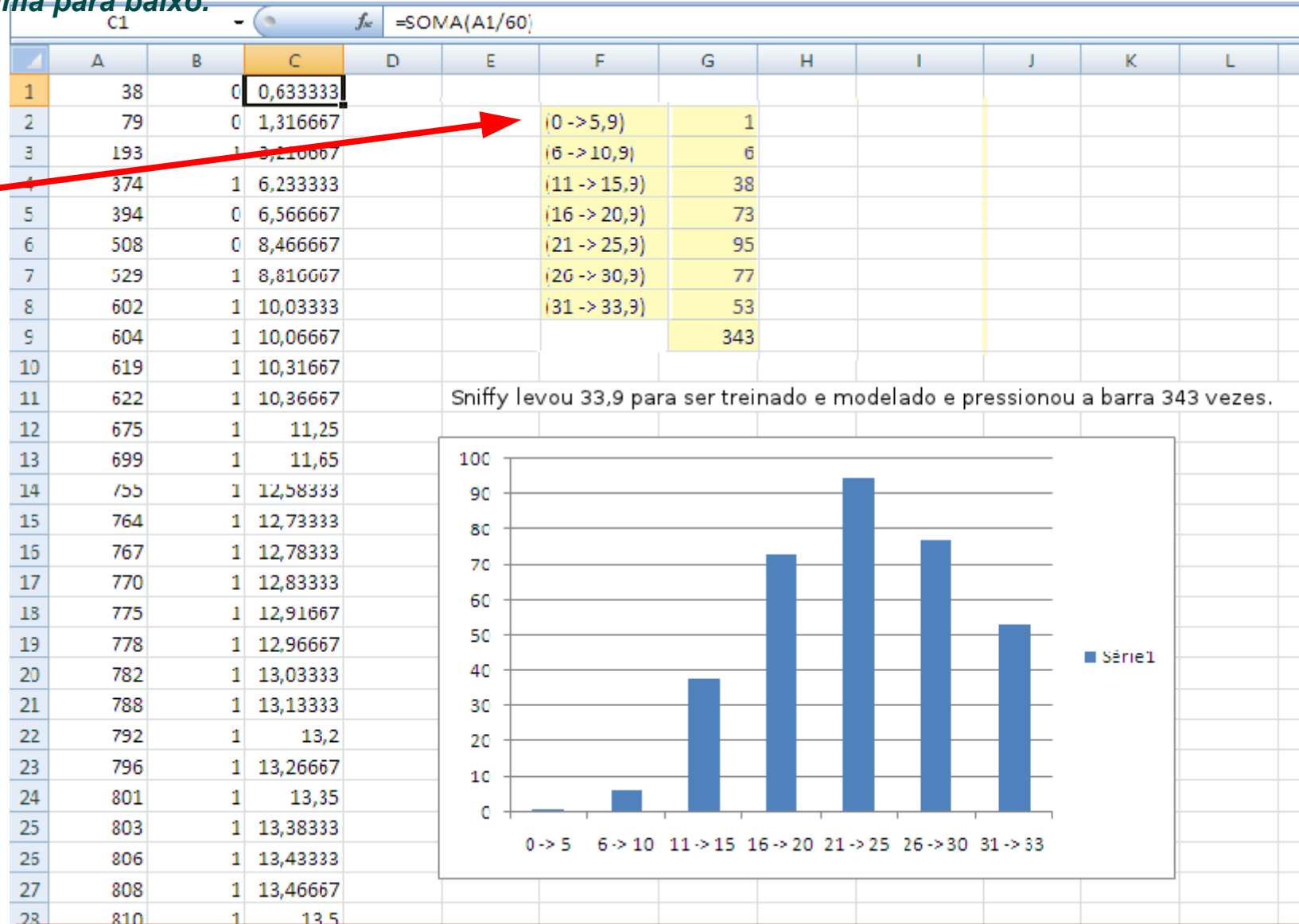
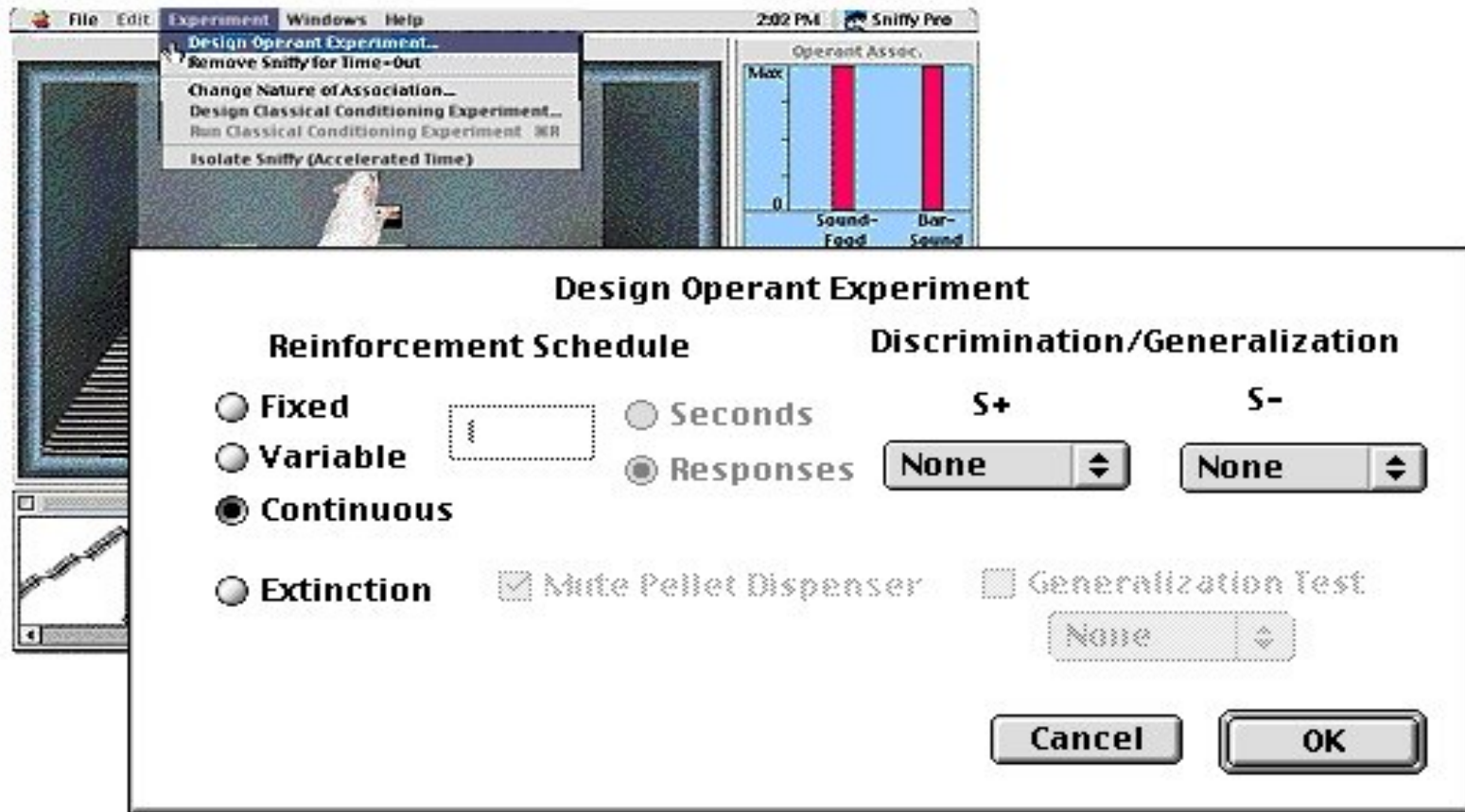
Extinction. What happens when reinforcement stops? To set up and run an extinction experiment, start with your trained Sniffy from the previous exercise (or OPEN the CRF Sniffy Data File in the Sample Files folder).

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
Sniffy will be pressing the bar frequently and eating after every press. When you choose the Design Operant Experiment command from the Experiment menu, you will see the dialog box shown above.

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

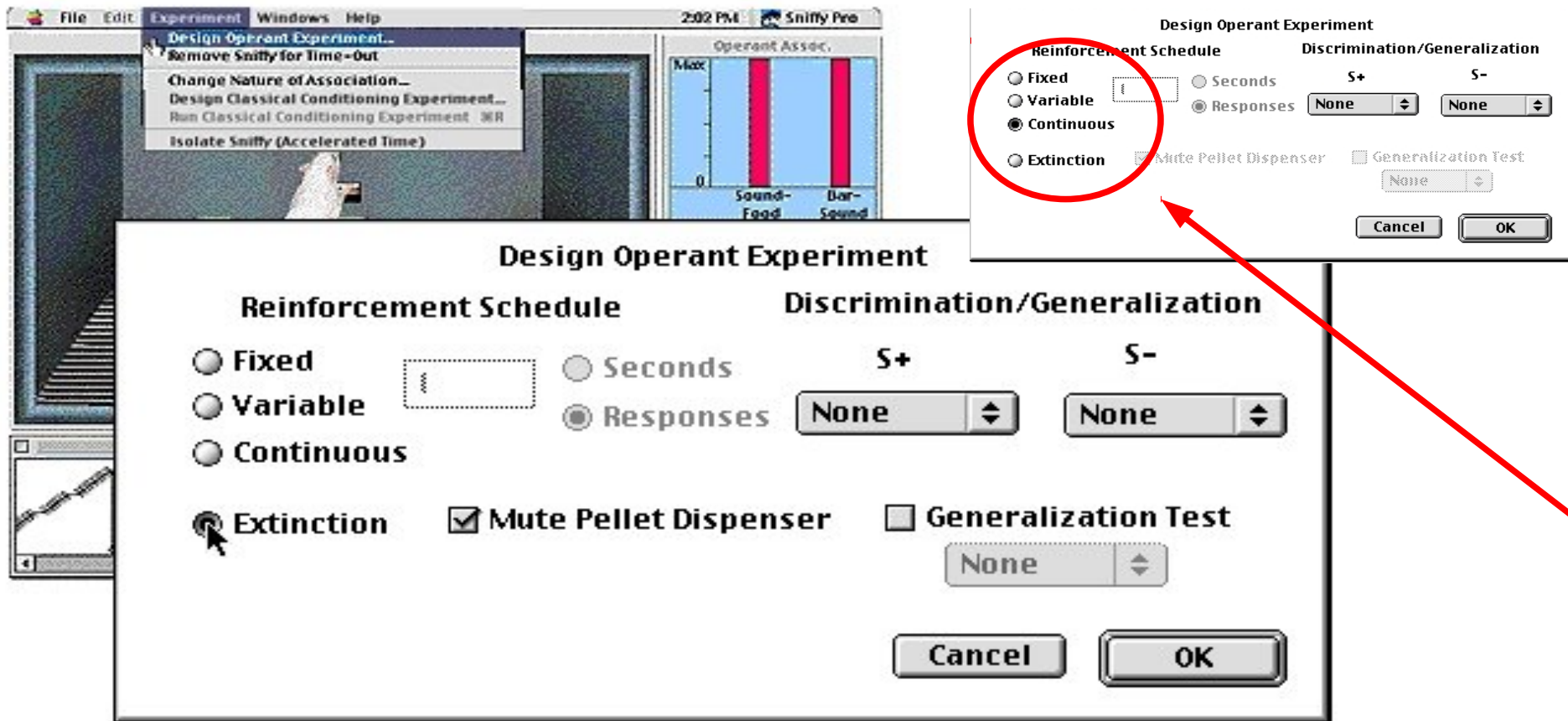
The button labeled "Continuous" is selected because every response has been reinforced; a procedure called continuous reinforcement. To select extinction, point the cursor at the button labeled "Extinction" and click your mouse button. Then click OK.

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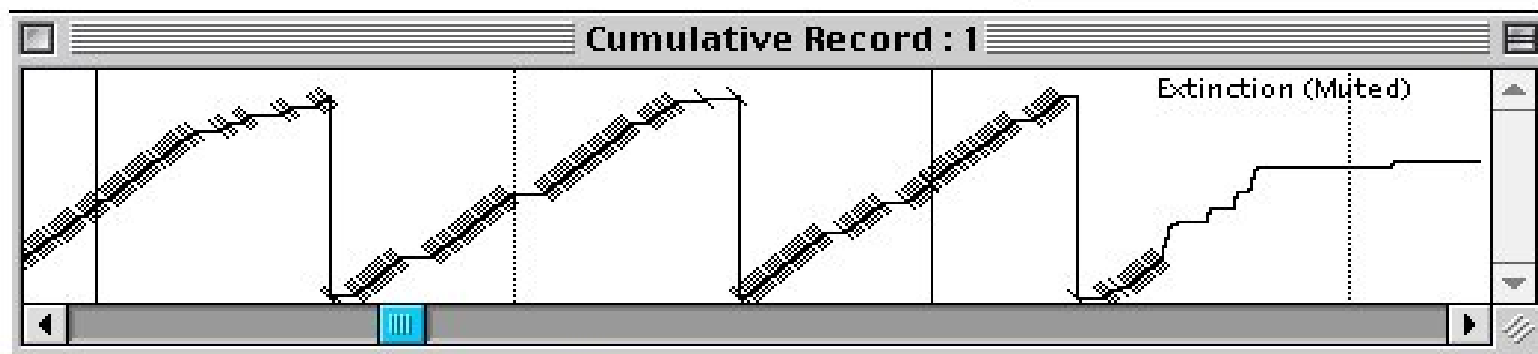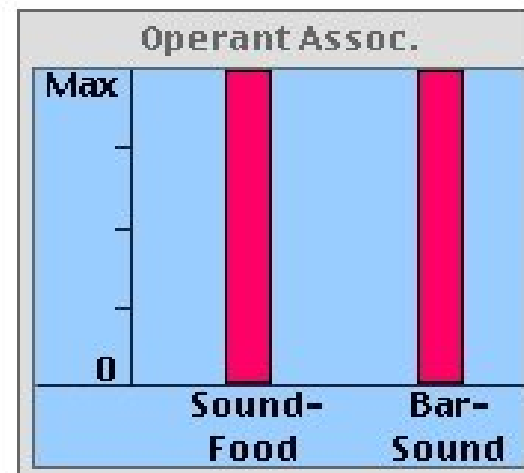
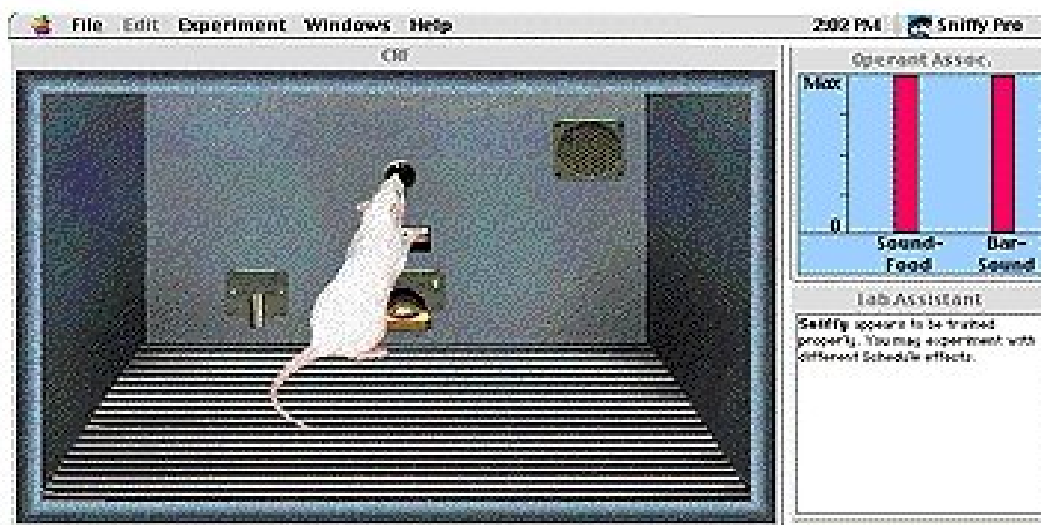
The program will begin running again, but Sniffy's bar presses will no longer be reinforced. Your cumulative record will look something like that shown above. Eventually the cumulative record will flatten out as Sniffy's rate of bar-pressing declines.

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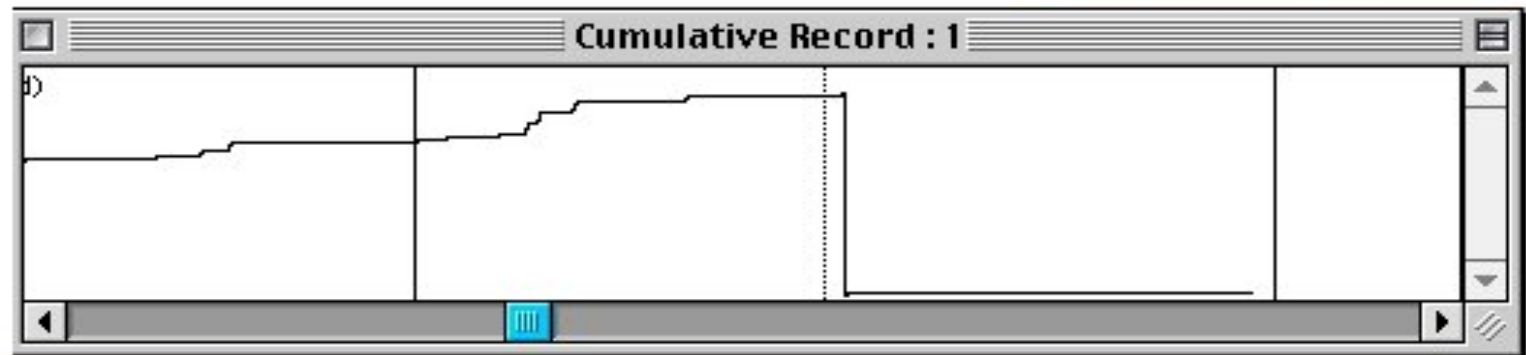
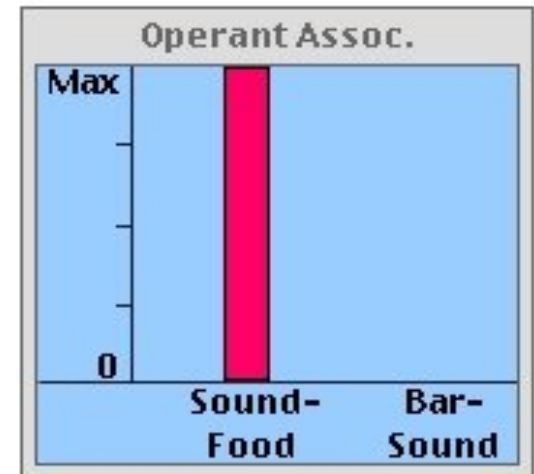
The extinction criterion is a 5-minute period during which Sniffy presses the bar no more than twice. At that point, save your Extinction file, and estimate the number of responses that Sniffy made and the time required to reach the extinction criterion.

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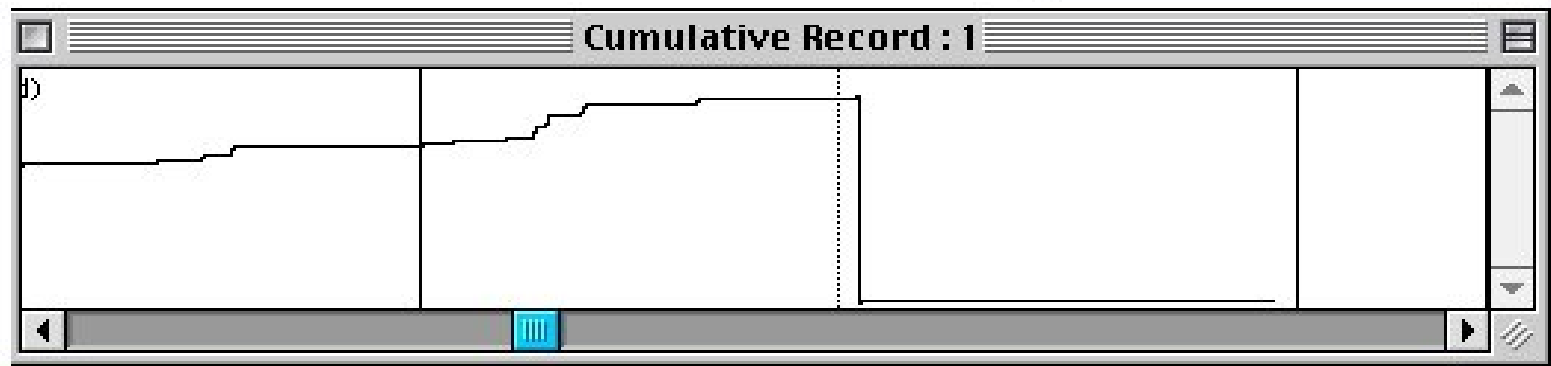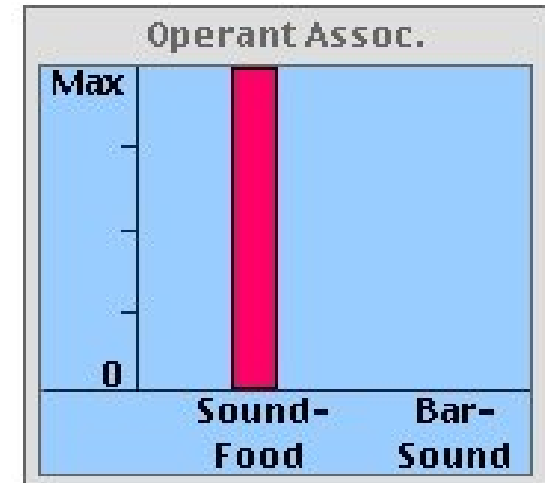
To measure the time required for extinction, you may have to scroll through the cumulative record and use the 5-minute lines as references. Watch as the cumulative record scrolls back in time.

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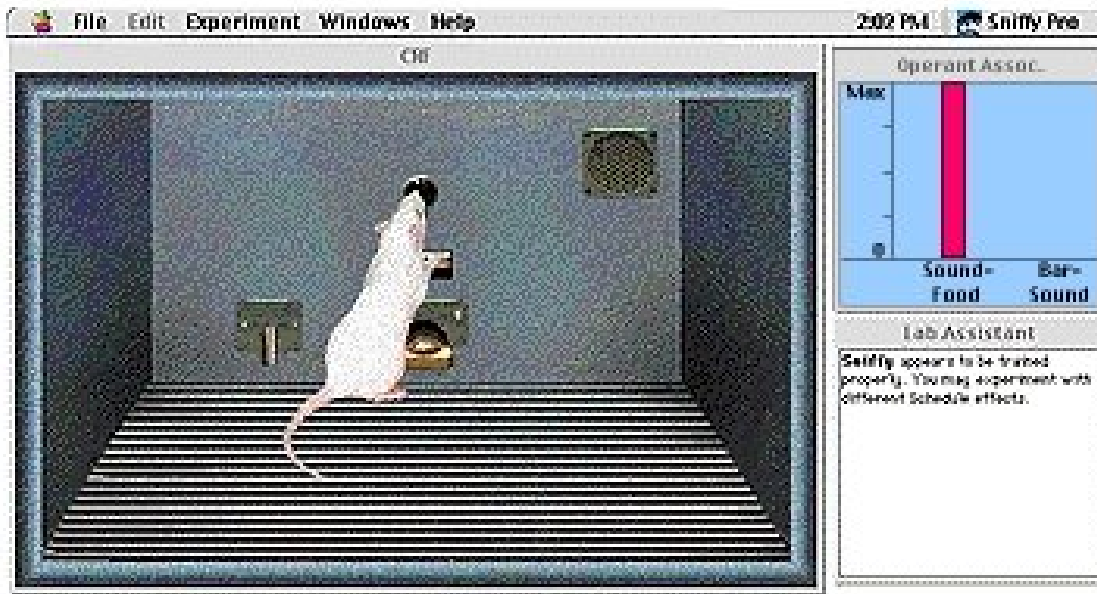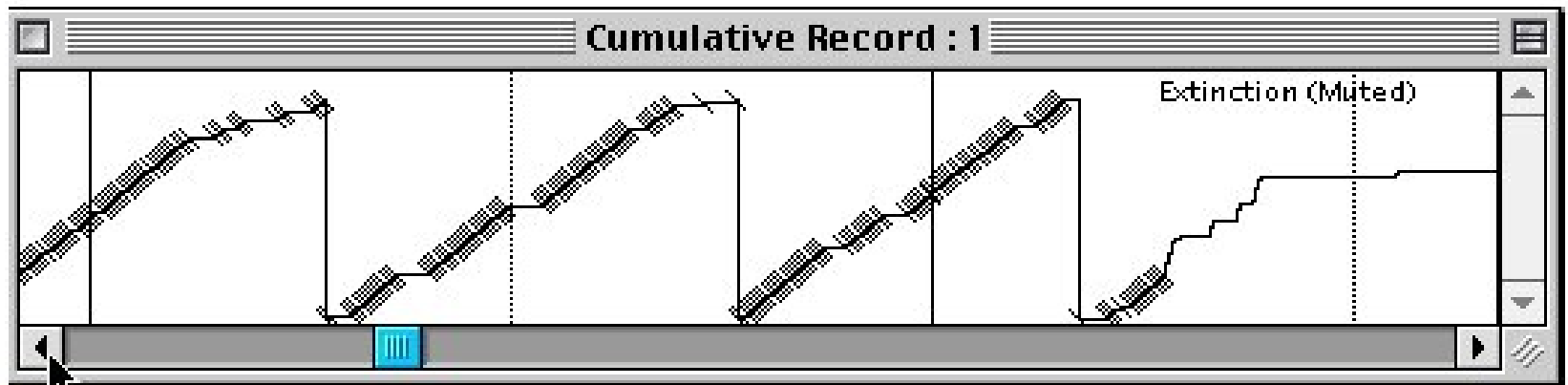
When estimating the time required for extinction, remember that alternating dotted and solid vertical lines mark off 5-minute periods. When estimating the number of bar presses, remember that Sniffy always makes 75 responses between two successive pen resets.

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
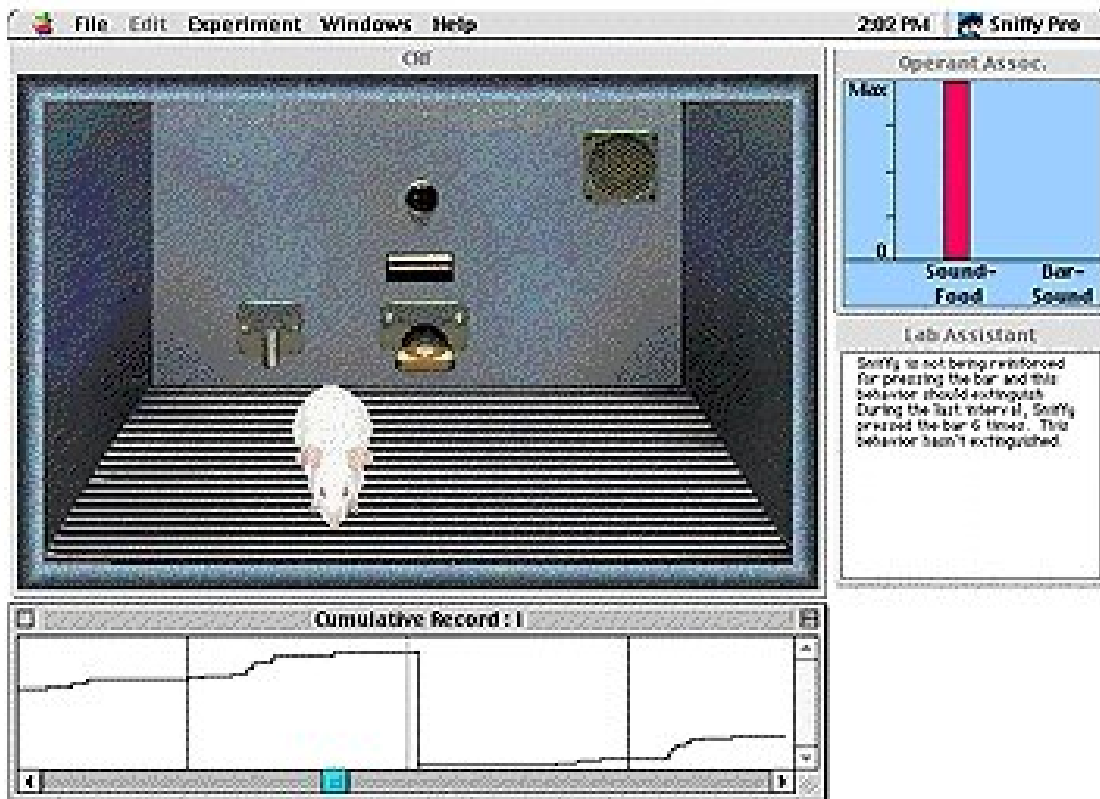
Spontaneous Recovery. A single extinction session will not permanently eliminate an operant response. If the rat is removed and returned 24 hours later, the operant response reappears. This rest-produced reappearance of an extinguished operant response is called spontaneous recovery.

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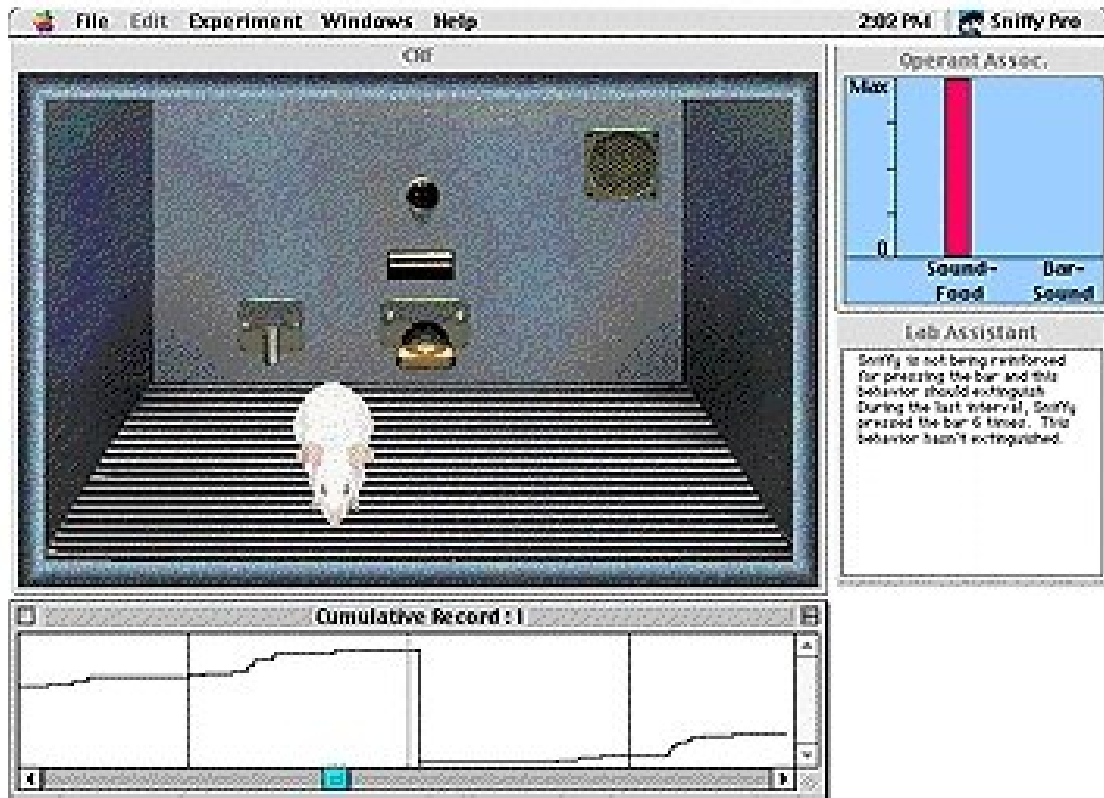
To simulate spontaneous recovery, start with a file in which Sniffy has been trained to press the bar and then extinguished. Then choose Remove Sniffy for Time-Out... command from the Experiment menu.

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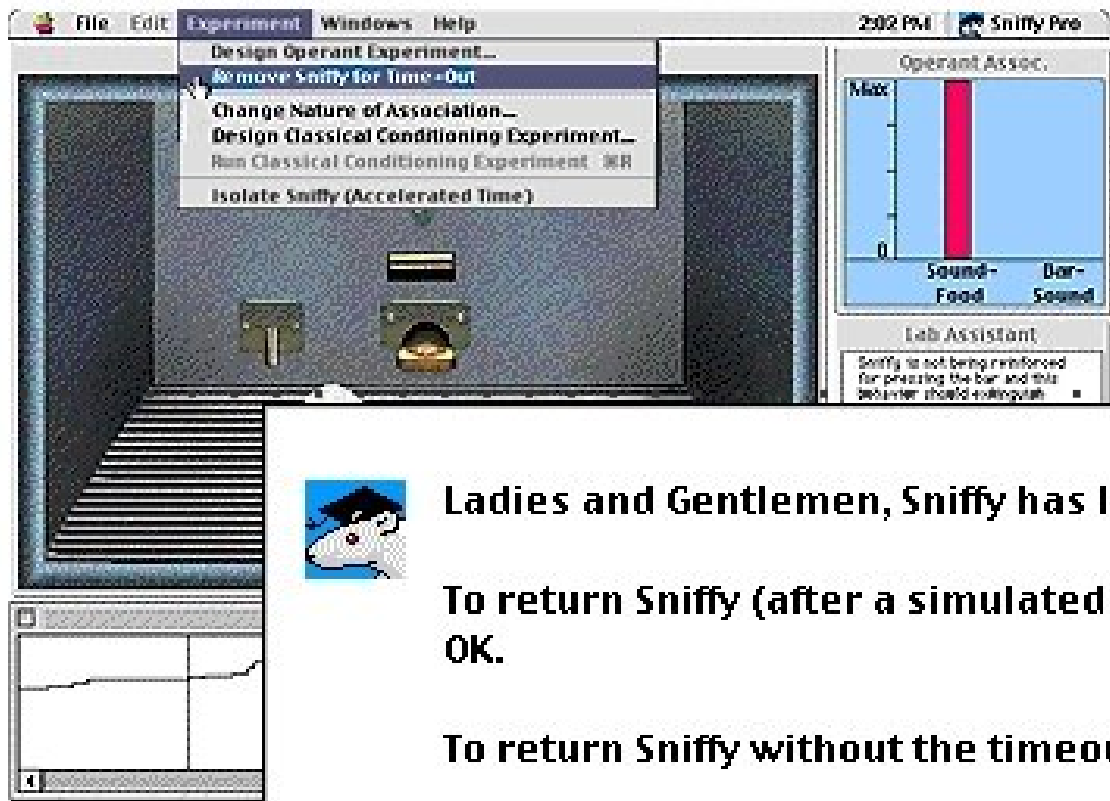Ladies and Gentlemen, Sniffy has left the chamber!

To return Sniffy (after a simulated time-out), click on OK.

To return Sniffy without the timeout, click on Cancel

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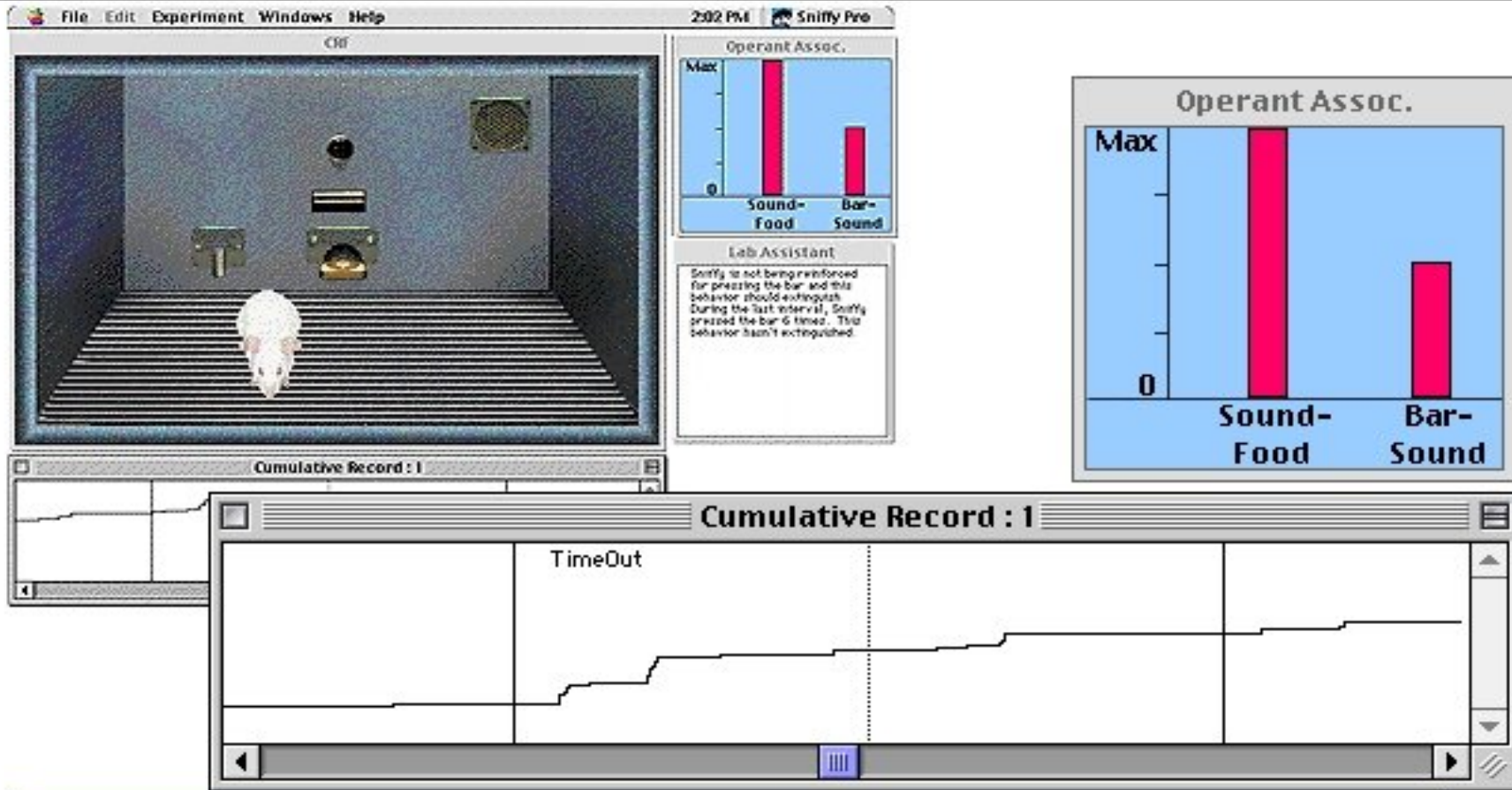
To simulate taking a rest, Sniffy will briefly disappear from view. When he reappears, he will temporarily recover bar-pressing activity. Immediately after the time-out, the cumulative record will resemble that shown next.

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
Let the program run until the extinction criterion is met again. Initially, the Operant Associations window shows the extinguished association is partly re-established. This is the "psychological" reason why Sniffy temporarily presses the bar again at the beginning of the second extinction session.

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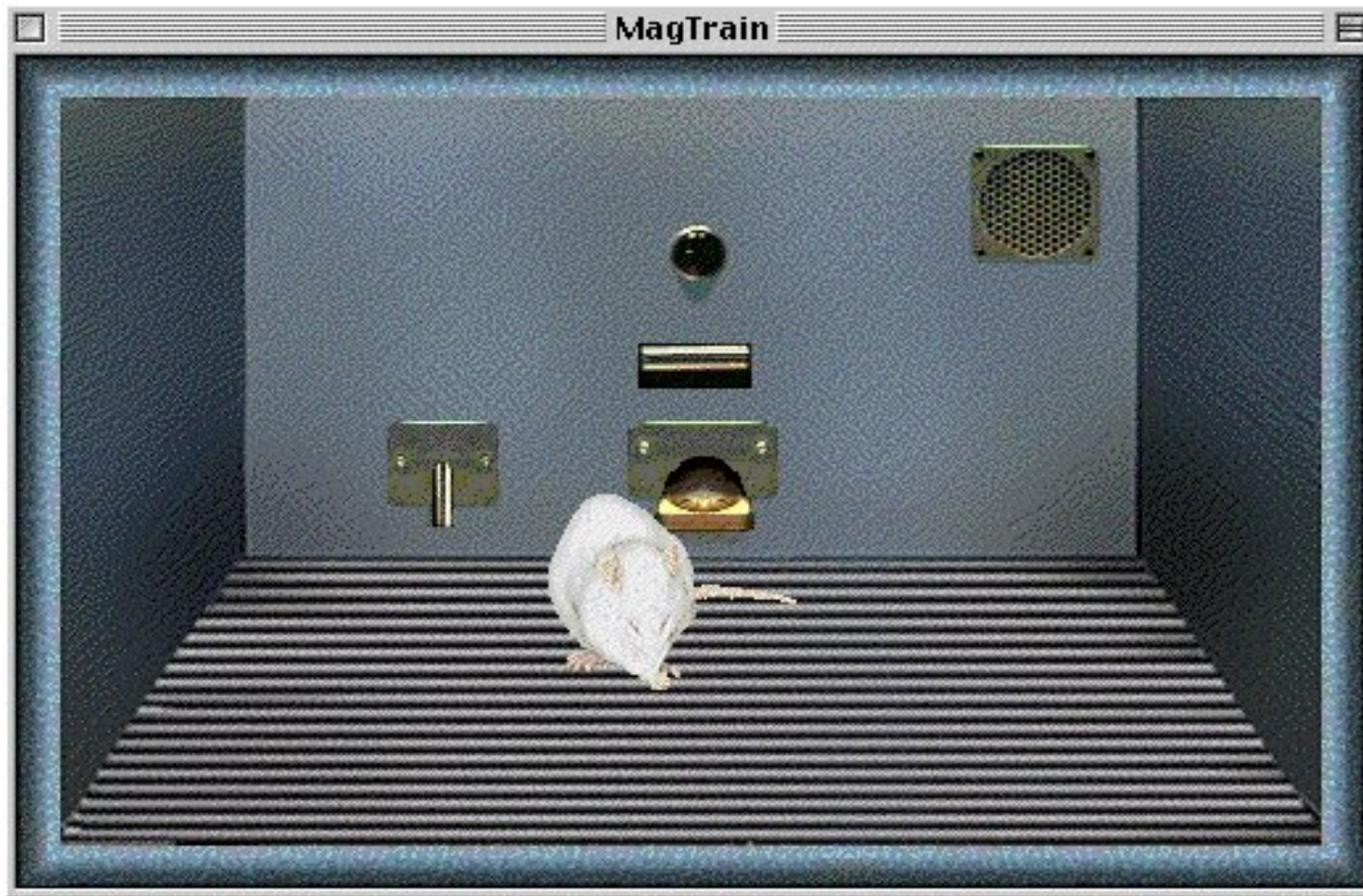Submitting your work. Save your file with an appropriate new name (e.g. SponRec) after the second extinction session. Your instructor may ask you to submit your file on a floppy disk or as an e-mail attachment.

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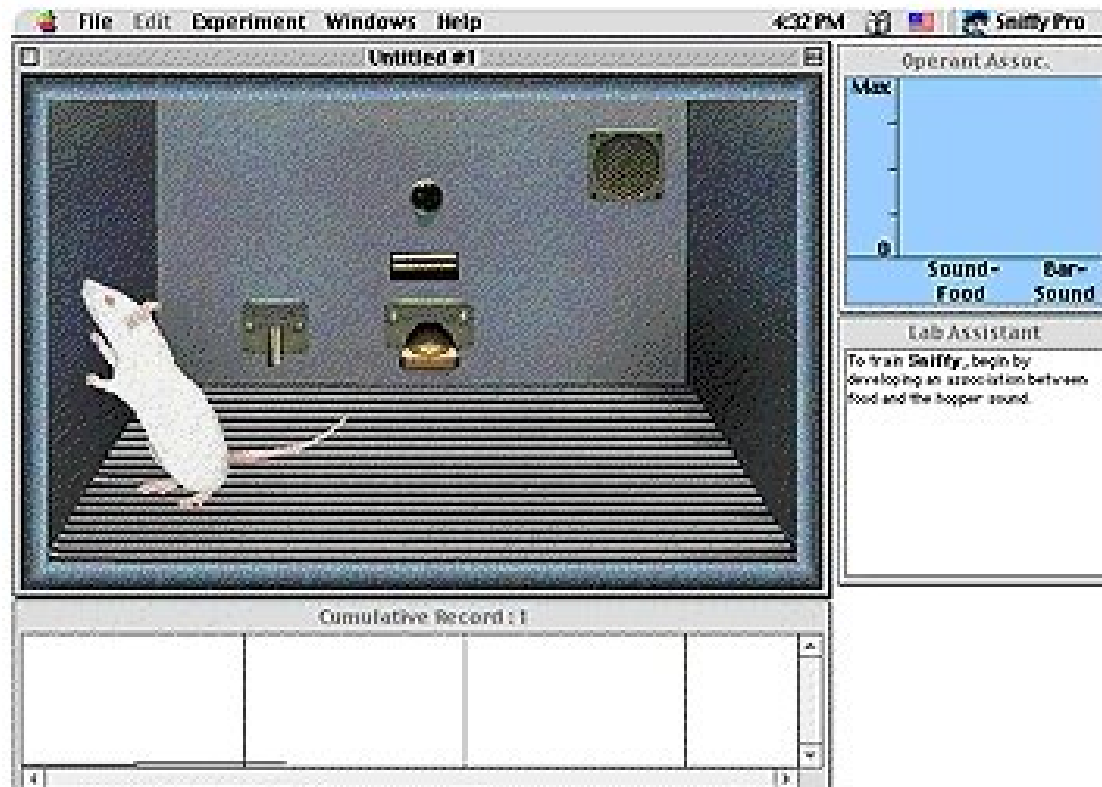
Thank you for exploring the
first Project. See Projects 2
and 3 for more advanced
material.

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